


U-Probe Grounding & Referencing Jumper Instructions


The U-Probe is shipped with one wire-attached jumper and one bare jumper. Use the wire-attached jumper for grounding metal components on or around the animal.

8- or 16-channel U-Probe:


1) To leave the tube on the U-Probe floating, use the following jumper configuration:

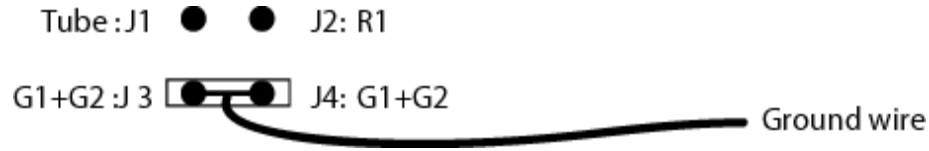

2) To ground the tube on the U-Probe, use the following jumper configuration:

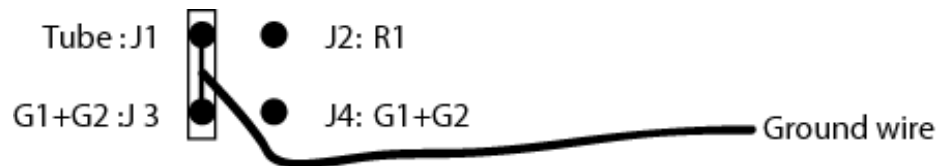


3) To attach the tube to the Reference (Ref 1) pin on the connector, use the following jumper configuration:

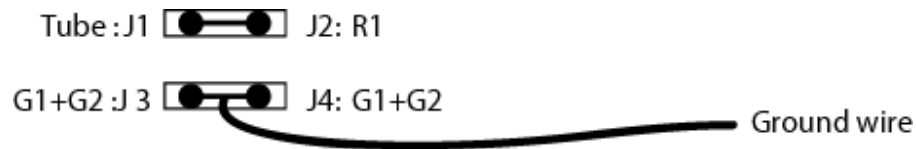


24-channel U-Probe:

1) To leave the tube on the U-Probe floating, use the following jumper configuration:

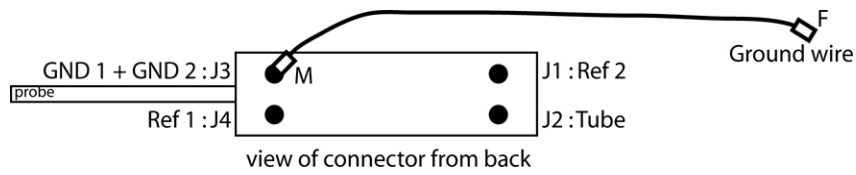

2) To ground the tube on the U-Probe, use the following jumper configuration:

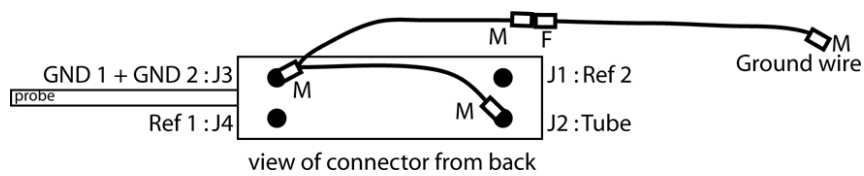

3) To attach the tube to the Reference (Ref 1) pin on the connector, use the following jumper configuration:

